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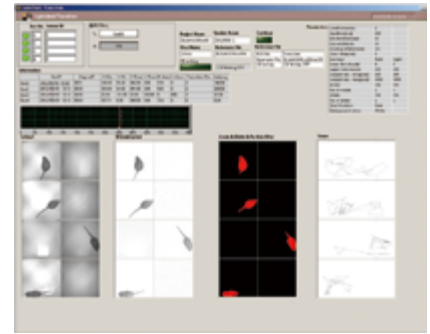
Light/dark transition test <emotionality>

Conditioned place preference test <drug dependency>

With added components, you can perform two types of experiments in this apparatus; light/dark transition test (2 compartments) or conditioned place preference test. In the light/dark transition test, one option is to customize the slide door connecting the two compartments to be automated and controlled by the software. Furthermore, the ceiling of each compartment has CCD camera in which images are analyzed by the software to measure the subject's movement within each compartment. In addition to number of entries and cumulative duration parameters, distance traveled within each box can also be measured. In conditioned place preference test, the light/dark compartments are recommended to be placed in sound-attenuating chambers.

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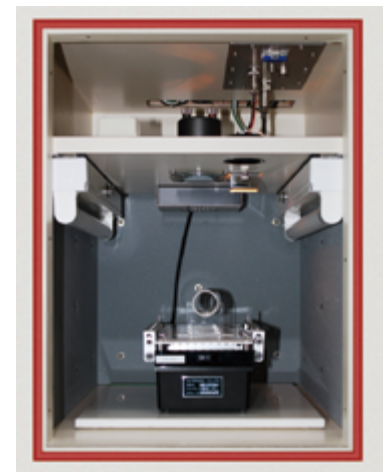
<センサーモーターゲーティング・短期的馴化・長期的馴化>

Startle response & Pre-pulse inhibition test

<sensorimotor gating / short-term habituation / long-term habituation>

※Multi chamber system (1 – 8 chamber)

On this system, you will put subject into cylindrical chamber and put on an acceleration sensor, then software quantify the startle response of subject. When you use chamber for this test, you need to put the chamber in sound attenuating box. You can easily set or modify sensitivity of acceleration sensor with attached gauge. By using software, you can manage schedule arbitrarily, save measured data, and check the data as waveform. Also, software saves the data as text format.

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