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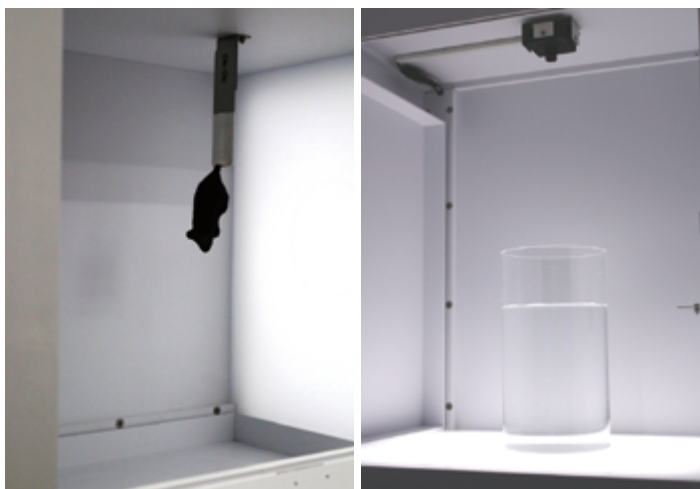
Porsolt forced swim & Tail suspension test
<emotionality ; depressant / antidepressant test>

※Multi chamber system (1 – 4 chamber)

Both Porsolt forced swim test and Tail suspension test, freezing can be automatically measured by image analysis. Images of subjects in the forced swim test are captured from directly above, while subjects in the tail suspension test are taken from the side. Lighting within the chamber illuminates the background of the subject, thus decreasing the amount of noise and allowing for a clear image analysis. Furthermore, the tail suspension test has been designed to prevent the subject from climbing up on its own tail during testing to provide more accurate behavioral measurements. マウスの強制水泳実験とテールサスペンション実験が可能なシステムです(ラットは強制水泳実験のみ)。動物の背景にLEDパネルライトを設置して撮影し、その動画をリアルタイムで画像解析することで、動物の無動時間を自動的に数値化します。1~4個体同時に実験可能です。

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Vogel type conflict test <emotionality / anxiolytic test>

※Multi chamber system (1 – 5 chamber)

On this system, software measures number that subject licks tube valve of a water bottle. Putting water-deprived subject into chamber and let it drink water at first, and every time subject licks valve certain times you set, electrical shock is provided through its mouth. You can set licking number to provide electrical shock and strength of shock. By using software, you can manage schedule and measure data automatically.

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