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Three-points Meter

Activity, food intake and water intake monitor

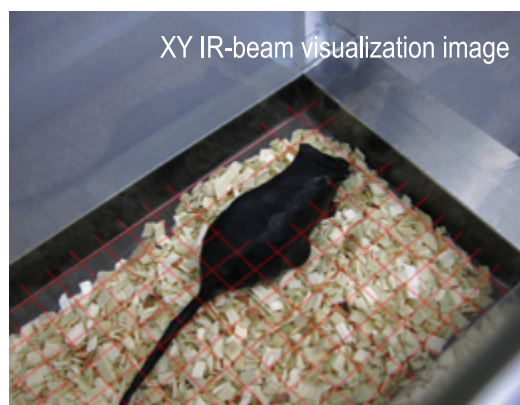
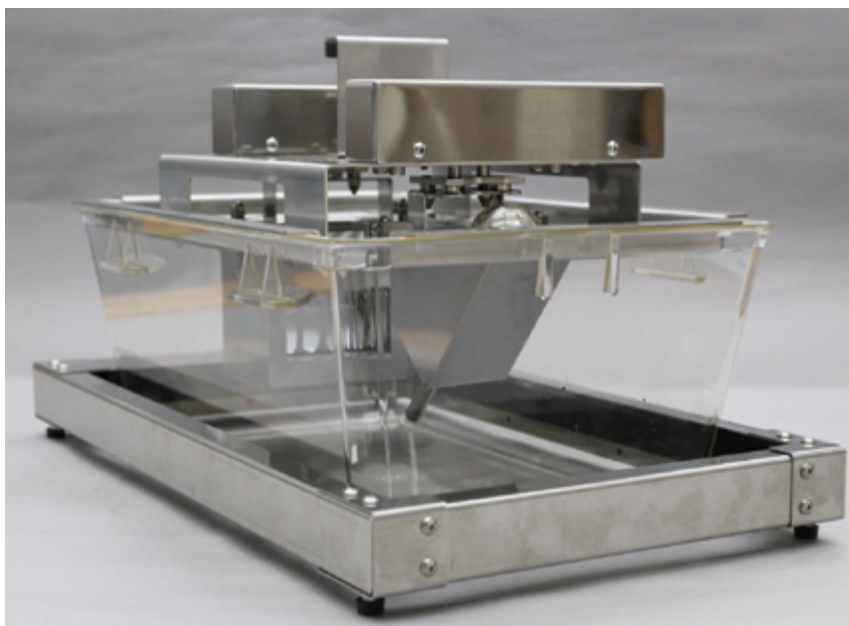
<amount of activity, food intake and water intake / biological rhythm >

※Multi chamber system (1 – 16 chamber)

Activity data acquisition software

Food and Water intake data acquisition software

This system can measure for 3 kinds of amounts: 1 ambulatory activity, 2 food intake and 3 water intake, at the same time within one cage. It is possible to measure at the same time such amount within 16 cages at the maximum (16 individuals) by means of a computer. For measuring activity, the position of the target is detected by interrupting the infrared beams positioned in the X and Y axis around the cage. This device detects the movement in position of the whole body of the target animal, without being misled by the movement of the tail and any other body. The position and movement of the target within the cage are automatically measured. The food box and the water bottle are placed under a precision scale and the software periodically measures the weight of the box.

Food intake and water intake are calculated by the difference between the measured weights. The apparatus is designed to eliminate dropped food, bedding material, and defecations to be added into the measurements. By using the software, time series data can automatically be saved as text files.

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